



Aliquot Preparation Kit



Figure 1: Aliquot Preparation Kit

The Aliquot Preparation Kit (Figure 1) contains the following components:

Component	Quantity
Silicon Oil	1 Bottle
Silicon Oil Stamp (standard \varnothing 6 mm)	1 Unit
Template for Silicon Oil Stamping	1 Unit
Aliquot Storage Trays	5 Units
Storage Case	1 Unit



Technical Specification

Please find attached the technical datasheet or contact us.

How to use the Aliquot Preparation Kit?

Step 1: Open the screws on both ends of the Aliquot Storage stack to reveal the Aliquot Storage Trays (Figure 2).



Figure 2: Aliquot storage

Step 2: Unbox the Template for Silicon Oil Stamping.

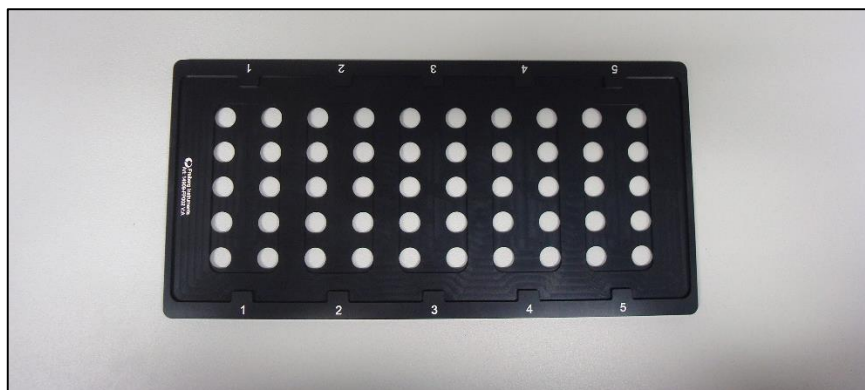


Figure 3: Template



Step 3: Place the suitable sample cups on the Aliquot Storage Trays for Silicon Oil Stamping and then place the Template for Silicon Oil Stamping on top of the Aliquot Storage Tray as shown in Figure 4.



Figure 4: Sample cups added to the Aliquot Storage Trays

Step 4: As shown in Figure 5, stamp up to 10 sample cups with Silicon Oil at a time.



Figure 5: Using the Silicon Oil Stamp

Step 5: Remove the Template and place the Aliquots on the sample discs.